

Ladd Research
83 Holly Court
Williston, VT 05495
Tel: (802) 658-4961
Fax: (802) 660-8859
Email: sales@laddresearch.com
Web: www.laddresearch.com

INSTRUCTION SHEET

SAMPLE HOLDER FOR THE HITACHI SEM 4800 (SERIES I OR II)

This H4800 is available in two versions:

Series I
Series II

The height requirement varies on the Series I and Series II

4800I - This is the most popular. It requires a height of 34mm to 36mm, measured from the underside of the "T" nut.

4800II - This requires a height of 29mm to 27mm.

The **LADD** holder is adjustable to fit the series I and II. By removing the 7mm spacer which is built into the holder.

To remove spacer: a) Remove the screw from the base which holds the spacer.

b) Then reassemble the holder without the spacer.