

Ladd Research
3 Ewing Place
Essex Junction, VT 05452
Tel: (802) 658-4961
Email: sales@laddresearch.com
Web: www.laddresearch.com



Bell Shine

Catalog Number: 30099

Instructions for Use

1. Spray 2 or 3 bursts of Bell Shine onto interior of bell jar.
2. Using lint-free cloth, thoroughly wipe inside of bell jar to spread a very thin coating of Bell Shine.
3. When ready to clean, spray bell jar with cleaner (e.g., Windex).
4. Wipe bell jar clean with dry paper towels.
5. Rinse with damp paper towels until the cleaner is removed.
6. Dry bell jar.
7. Recoat with Bell Shine before reuse.